HANDLEIDING - MODE D'EMPLOI - MANUAL

UCI120 (754750221)

UCI175 (754750222)

Ultrasoonreiniger Nettoyeur à ultrasons Ultrasonic cleaner

- NL** P.02 Gelieve te lezen en voor later gebruik bewaren
- FR** P.07 Veuillez lire et conserver pour consultation ultérieure
- EN** P.12 Please read and keep for future reference

1 Veiligheidsvoorschriften

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1 Safety instructions

- Please read this instruction manual carefully before use and keep it for reference.
- Ensure correct power and switch connection before starting the equipment.
- Be careful the control panel will be corroded under the condition of organic solution, strong acid or strong base.
- Do not use the product under the following conditions:
 - Where temperature fluctuate strongly,
 - Where it has high humidity and especially dew,
 - Where vibration or shock is strong,
 - Where corrosive gas or dust exist,
 - Where water, oil or chemicals splash,
 - Where is easy to cause explosion.
- Don't let the liquid splash into the electronic parts, especially transducers.
- Don't contact the bottom directly (vibration plate), or may damage the transducers.
- Keep the machine in dry and cool environment.
- The machine power supply must be grounded.
- Wash the tank if there is much deposit in the tank
- Don't move the machine when there is liquid inside tank to avoid splash.


2 Technical data

Model	UCI120	UCI175
Voltage	3x400 V	220 V
Tank capacity	120 litre	175 litre
Ultrasonic power	1800 W	2400 W
Heating power	3000 W	4500 W
Liquid temperature setting	20 °C - 80 °C	20 °C - 80 °C
Tank dimensions	600 x 500 x 400 mm	700 x 500 x 500 mm
Timer	1 - 99 minutes	1 - 99 minutes

3 Cautions before starting the machine

- Place all buttons on OFF.
- Fill the cleaning liquid tank before switching on the machine. Do not empty the tank when machine is working.
- Start the ultrasonic cleaning when the liquid has reached the set temperature.
- Don't start the heating if there is no liquid or the liquid level is lower than the work level, or heating pad will be damaged.
- Put the lid on the tank to reduce the noise and to avoid liquid loss and evaporation.

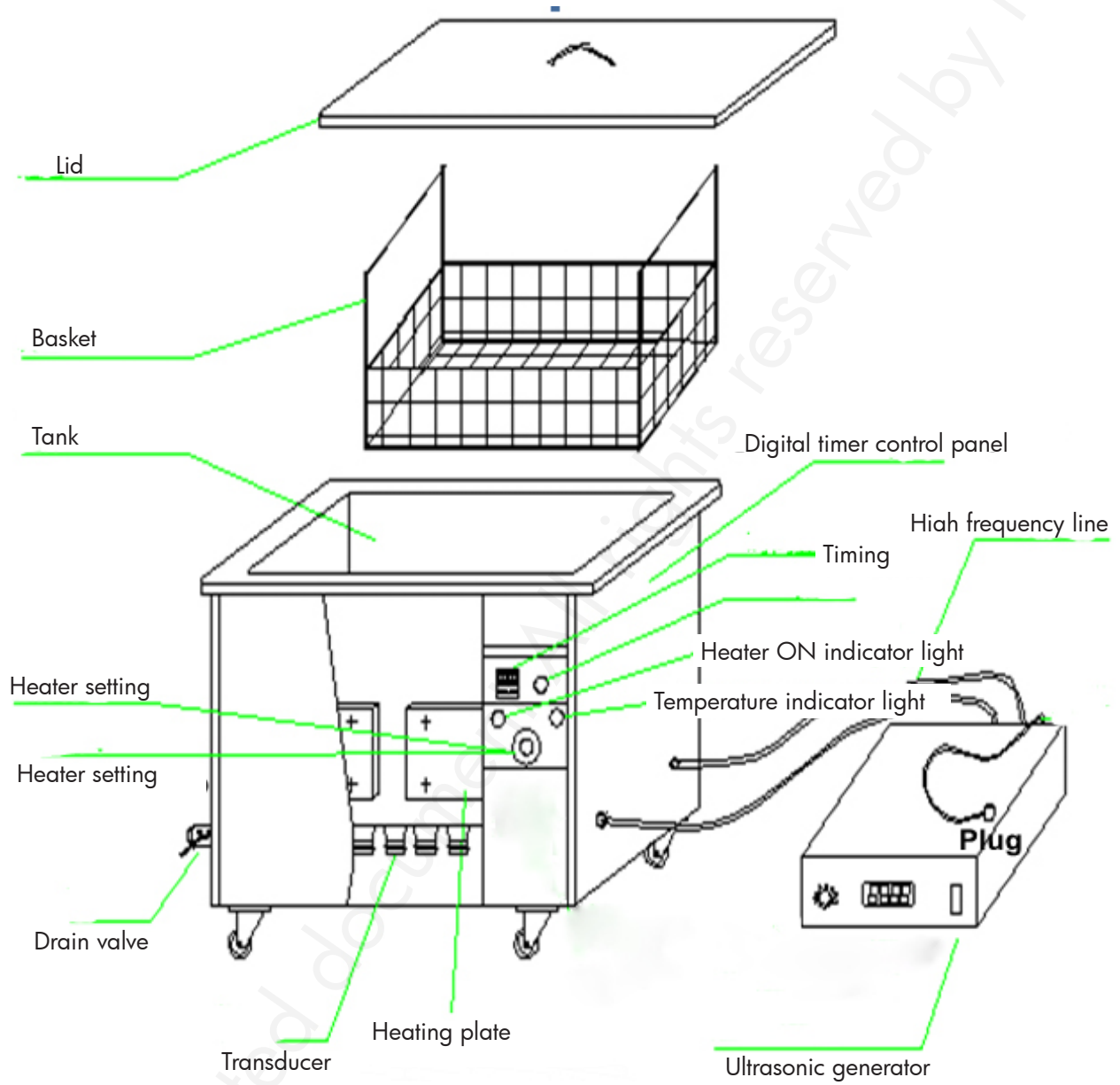4 Structure diagram

Generator separated control

5 Operation

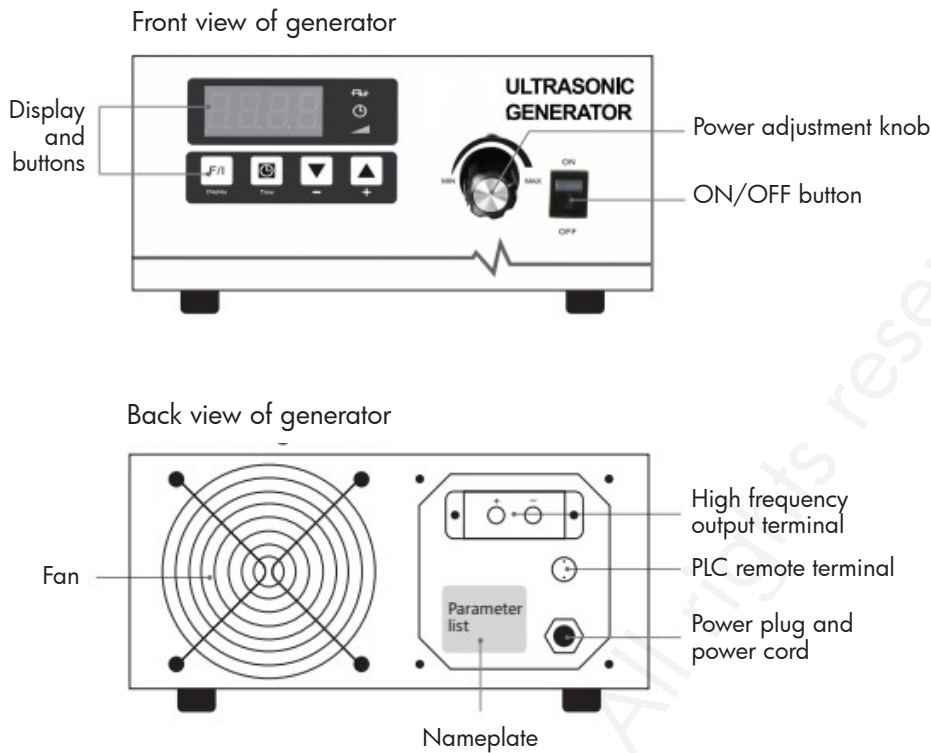
5.1 Control panel

Control panel

- **ON/OFF button**
To start or stop ultrasonic.
- **Temperature adjustment knob**
Turn the knob clockwise to increase temperature, turn the knob counterclockwise to decrease temperature.
- **Heater ON**
Heater working indicator. It turns red when heating system is working.
- **Temperature reach**
Heater indicator, it turns green when temperature reaches the setting value, heating system will stop working.
- **Timer display**
Shows ultrasonic working time. Maximum working time is 99 hours. Time can be set by hour or by minute.

5.2 Ultrasonic generator

5.3 Operating instructions

1. Place the machine on a flat floor.
2. Connect the power cable correctly.
3. Connect the drain pipe with factory drainage system.
4. Put the cleaning liquid into the tank.
5. Immerse the workpiece into the tank basket.
6. Set the temperature based on requirements and turn on the heater switch.
7. When the setting temperature is reached, adjust the ultrasonic power to Min., then start ultrasonic and adjust ultrasonic power to requirement after normal vibration.

6 Maintenance

- Use high-pressure blast to clean dust of the control system regularly.
- Clean the tank once a week at least.

7 Trouble shooting

Problems	Possible causes	Solutions
No ultrasonic.	Machine not connected.	Check and connect the machine.
	Fuse broken.	Check power supply and fuse. Replace it if necessary.
	Cable short circuit.	Replace power cord.
	Transducer short circuit.	Contact the technical service of your dealer.
	Other reasons.	Contact the technical service of your dealer.
Unsatisfactory cleaning.	Ultrasonic cleaning not strong enough.	Adjust ultrasonic button.
	Liquid level too high or too low.	Adjust liquid to the best level.
	Temperature too high or too low.	Adjust temperature to the most appropriate level. Suggestion: 50-60 °C.
	Cleaning liquid not suitable.	Stop and switch off the power supply, replace liquid after the previous liquid cooled down.
	Other reasons.	Contact the technical service of your dealer.
No heating.	Bad connection of heating power switch.	Check heating plug to correct connection.
	Fuse broken.	Check out the line with multimeter: if OK and resistance value is few hundreds then replace the fuse. If not OK, it's short circuit, replace heater.
	Other reasons.	Contact the technical service of your dealer.
Temperature control failure.	Thermostat loosen.	Fasten thermostat header.
	Thermostat tube broken.	Replace thermostat.
	Other reasons.	Contact the technical service of your dealer.
Timer control failure.	Timer knob out of control.	Loosen or tighten the screw.
	Timer failure.	Replace timer or digital panel.
	Other reasons.	Contact the technical service of your dealer.
Electric leakage.	Electrical outlet not grounded.	Ensure proper grounding of electrical outlet.
	Machine not grounded.	Ensure proper grounding of the machine.
Other problems.		Contact the technical service of your dealer.

- NL** 8 **EG conformiteitsverklaring**
FR 8 **Déclaration de conformité CE**
EN 8 **EC declaration of conformity**

Fabrikant/Invoerder
 Fabricant/Importateur
 Manufacturer/Retailer

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Meets the provisions of the aforementioned directive, including, any amendments valid at the time of this statement.

Mouscron, 20/09/2021

Bart Vynckier, Director
 Vynckier Tools sa